

Determination of Nitrogen in Flour

LECO Corporation; Saint Joseph, Michigan USA

Instrument: TruMac[®] N

Introduction

Flour is a fine particle powder created by milling or grinding a dry grain. The most common varieties of flour are made from wheat, although any grain can be used to make flour. Flour is typically used to make dough for a variety of bread products. The protein content in the flour is one of the primary constituents that determines the best use for the flour, with lower-protein flour (~8%) typically being used for cakes and pastries, mid-range protein flours (~10%) being categorized as all-purpose, and higher-protein flours (~12%) being referred to as bread flour.

The accurate and precise determination of protein not only plays a role in the characterization of nutritional or dietary value in flours, but is also the key to determining the category or quality of the flour. Protein in flour and other food products is most commonly calculated using the measured nitrogen in the sample and a protein factor multiplier (protein factors vary according to the sample matrix).

The LECO TruMac N is a macro combustion nitrogen/protein determinator that utilizes a pure oxygen environment in a ceramic horizontal furnace and large ceramic boats for the macro sample combustion process. A thermoelectric cooler removes the moisture in the combustion gas without the use of chemical reagents. A 3 or 10 cc volume of combustion gas is taken using a combustion gas collection and handling system. The combustion gas collection and handling system achieves a low cost-per-analysis by reducing the amount of chemical reagents used for scrubbing and converting the nitrogen oxide combustion gas to nitrogen. A thermal conductivity (TC) cell is used for the detection of nitrogen in the combustion gas.

Sample Preparation

Samples must be of uniform consistency to produce suitable results.

Accessories 528-203 Crucibles

Calibration Samples

502-092 EDTA, 502-642 Phenylalanine,
501-050 Nicotinic Acid

Analysis Parameters*

| | |
|-----------------------|-----------|
| Furnace Temperature | 1100°C |
| TE Cooler Temperature | 5°C |
| Dehydration Time | 0 seconds |
| Purge Cycles | 2 seconds |



Instrument Model and Configuration

Thermal conductivity detectors work by detecting changes in the thermal conductivity of the analytical gas compared to the constant thermal conductivity of the reference gas. The greater the difference between the thermal conductivity of the carrier gas and the analyte gas, the greater the sensitivity of the detector. The TruMac is available in models that support either the use of helium or argon as the instrument's carrier gas for the thermal conductivity cell.

When used as a carrier gas, helium provides the highest sensitivity, providing the best performance at the lower end of the nitrogen range. Helium models also offer the additional advantage of replacing the 10 cc aliquot loop with a 3 cc loop within the instrument's gas collection and handling system. The 10 cc aliquot loop optimizes the instrument for the lowest nitrogen range and best precision. The 3 cc aliquot loop extends reagent life expectancy by approximately three fold compared to the 10 cc aliquot loop, while providing the lowest cost per analysis with minimal impact on practical application performance (see Typical Results section).

Due to the recent history of low supply and general availability issues for helium gas, the argon model was developed to utilize argon as a carrier gas. Since the thermal conductivity difference between argon and nitrogen is not as great as the thermal conductivity difference between helium and nitrogen, the detector is inherently less sensitive with argon as a carrier gas. The argon model has a similar practical application performance compared to the helium model, operating with equivalent instrument and method configurations (see Typical Results section).

Note: Changing carrier gas and aliquot loop size requires hardware changes within the instrument.

Element Parameters

| | Helium | Argon |
|-----------------------|----------------|------------|
| | 10 cc and 3 cc | 10 cc |
| Baseline Delay Time | 6 seconds | 6 seconds |
| Minimum Analysis Time | 35 seconds | 55 seconds |
| Endline Time | 2 seconds | 2 seconds |
| Conversion Factor | 1.00 | 1.00 |
| Significant Digits | 5 | 5 |
| TC Baseline Time | 10 seconds | 10 seconds |

LECO[®]

Organic Application Note

Burn Profile

| Burn Cycle | Lance Flow | Purge Flow | Time (seconds) |
|------------|------------|------------|----------------|
| 1 | Off | On | 5 seconds |
| 2 | On | On | 35 seconds |
| 3 | On | Off | END |

Ballast Parameters

Ballast

| | |
|--------------------|-------------|
| Equilibrate Time | 30 seconds |
| Not Filled Timeout | 300 seconds |

Aliquot Loop

| | |
|---------------------------|-----------|
| Equilibrate Pressure Time | 4 seconds |
| High Precision | Yes |
| High Speed | No |

*Refer to TruMac Operator's Instruction Manual for Method Parameter definitions.

Procedure

1. Prepare instrument for operation as outlined in the operator's instruction manual.
2. Condition the system by analyzing 3 to 5 blanks (crucible is not required).
3. Determine blank.
 - a. Enter 1.0000 g mass into Sample Login (F3) using Blank as the sample name.
 - b. Place a 528-203 Crucible to the appropriate position of the autoloader.
 - c. Repeat steps 3a through 3b a minimum of three times.
 - d. Initiate the analysis sequence (F5).
 - e. Set the blank following the procedure outlined in the operator's instruction manual.

4. Calibrate.
 - a. Weigh ~0.75 g of EDTA calibration sample into a 528-203 Crucible, enter mass and sample identification into Sample Login (F3).
 - b. Transfer crucible to the appropriate position of the autoloader.
 - c. Repeat steps 4a through 4b a minimum of three times.
 - d. Initiate the analysis sequence (F5).
 - e. Calibrate the instrument following the procedure outlined in the operator's instruction manual. Use single standard calibration.

Note: Multi-point (fractional weight or multiple calibration samples) may be used to calibrate if desired. Research has shown that a properly functioning TruMac can be calibrated using several replicates of a single mass range (nominal 0.75 g) of EDTA utilizing a single standard calibration. This is a cost effective and simple process. The calibration can be verified by analyzing different compounds such as nicotinic acid (0.25 to 0.5 g) and/or phenylalanine (0.5 to 0.75 g).

5. Analyze Samples.
 - a. Weigh ~1 g of flour sample into a 528-203 Crucible; enter mass and sample identification into Sample Login (F3).
 - b. Transfer crucible to the appropriate position of the autoloader.
 - c. Repeat steps 5a through 5b for each sample to be analyzed.
 - d. Initiate the analysis sequence (F5).

Note: If soot (carbon black) is noticed in the primary filter (steel wool filter), reduce sample mass to prevent soot build-up in this filter. Soot can be produced when larger masses of some sample types are analyzed.

Typical Results

(based on single standard calibration using 0.75 g EDTA)

| | 3 cc Helium | | 10 cc Helium | | 10 cc Argon | |
|---------------------|-------------------|------|-------------------|------|-------------------|------|
| | Mass(g) | % N | Mass(g) | % N | Mass(g) | % N |
| LECO 501-563 | 1.0194 | 1.21 | 1.0070 | 1.21 | 1.0392 | 1.20 |
| Lot: 1014 | 1.0272 | 1.20 | 1.0069 | 1.20 | 1.0140 | 1.20 |
| 1.21±0.02 | 1.0221 | 1.20 | 1.0104 | 1.20 | 1.0401 | 1.20 |
| Corn Flour | 1.0490 | 1.20 | 1.0630 | 1.20 | 1.0289 | 1.20 |
| | 1.0275 | 1.20 | 1.0708 | 1.20 | 1.0091 | 1.20 |
| | Avg = 1.20 | | Avg = 1.20 | | Avg = 1.20 | |
| | s = 0.004 | | s = 0.004 | | s = 0.003 | |
| LECO 502-274 | 1.0658 | 2.74 | 1.0176 | 2.75 | 1.0038 | 2.74 |
| Lot: 1011 | 1.0209 | 2.74 | 1.0069 | 2.75 | 1.0389 | 2.75 |
| 2.75±0.02 | 1.0994 | 2.74 | 1.0107 | 2.76 | 1.0123 | 2.75 |
| Wheat Flour | 1.0157 | 2.75 | 1.0265 | 2.74 | 1.0135 | 2.74 |
| | 1.0091 | 2.75 | 1.0115 | 2.74 | 1.0377 | 2.74 |
| | Avg = 2.74 | | Avg = 2.75 | | Avg = 2.74 | |
| | s = 0.005 | | s = 0.008 | | s = 0.005 | |
| LECO 502-275 | 1.0044 | 1.73 | 1.0108 | 1.73 | 1.0118 | 1.72 |
| Lot: 1007 | 1.0080 | 1.72 | 1.0042 | 1.74 | 1.0312 | 1.72 |
| 1.74±0.06 | 1.0217 | 1.73 | 1.0194 | 1.74 | 1.0209 | 1.72 |
| Rye Flour | 1.0429 | 1.73 | 1.0421 | 1.73 | 1.0562 | 1.72 |
| | 1.0314 | 1.72 | 1.0114 | 1.73 | 1.0241 | 1.73 |
| | Avg = 1.73 | | Avg = 1.73 | | Avg = 1.72 | |
| | s = 0.005 | | s = 0.005 | | s = 0.004 | |
| LECO 502-278 | 1.0629 | 1.17 | 1.0465 | 1.17 | 1.0172 | 1.17 |
| Lot: 1013 | 1.0398 | 1.17 | 1.0182 | 1.17 | 1.0896 | 1.17 |
| 1.17±0.05 | 1.0384 | 1.17 | 1.0323 | 1.17 | 1.0170 | 1.17 |
| Rice Flour | 1.0560 | 1.17 | 1.0760 | 1.17 | 1.0689 | 1.17 |
| | 1.0220 | 1.17 | 1.0580 | 1.17 | 1.0371 | 1.17 |
| | Avg = 1.17 | | Avg = 1.17 | | Avg = 1.17 | |
| | s = 0.002 | | s = 0.001 | | s = 0.001 | |

TruMac is a registered trademark of LECO Corporation.

LECO Corporation | 3000 Lakeview Avenue | St. Joseph, MI 49085 | Phone: 800-292-6141 | Fax: 269-982-8977
 info@leco.com • www.leco.com | ISO-9001:2008 HQ-Q-994 | LECO is a registered trademark of LECO Corporation.