


LECO introduces the new PAX-it 2, featuring substantial improvements to the software layout and PAXCam Camera interface.

The new ribbon view ushers in a modernization that makes the system easier to use, and will help keep the user more organized. The context-sensitive nature of the software greatly declutters your workspace by hiding icons that are not in use.

Changes were also made to the browsing structure of the PAX-it image collection. Images and data can now be grouped more easily using the new tree view. At a glance, this tree view clearly shows your location in your project folders ,and allows for effortless navigation.

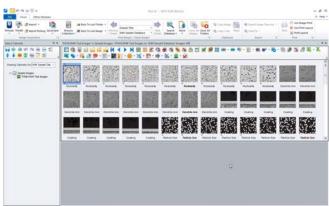

Select Cabinets 📮 🗙	'PAXit/IA44 Test	Images' is Samp	ole Images->PAX	t/IA44 Test Imag	es on 'IA44 Sam	ple Database' Ima	ages: 149							ą×
	📘 🐑 🖥 WK Wy - E	🧊 💼 🕼 • 🥷 🚇 1	i 🛶 성 ┥	📕 🖌 🖣	🌐 💋 🛍 (- 👧 -)))	🔕 🛸 🐜 - 🏡 - 🌨) 🛍 🏝 🗾 - 🖌	í 🗹 😰 🔡	🖶 - 🖶 i	🧚 • 📄 • 🔮	u @ - 🌄	- 🛍 🛃 🍒	/ 🚑 🖾 - X	¢ .
🕇 🖊 🔜 🖺 😔 🗾 圮 🖕	1													•
Viewing Cabinets On: Hard Drive														=
⊡ - [] {Untitled}	Nodularity	Nodularity	Nodularity	Nodularity	Nodularity	Nodularity	Nodularity	Nodularity	Nodularity	Nodularity	Nodularity	Dendrite Arm	Dendrite Arm	
Calibration Images														
Company A	Constant States								7 12 19 19 19 19			RUSECICE		
Lot #2	Dendrite Arm	Dendrite Arm	Dendrite Arm	Dendrite Arm	Dendrite Arm	Dendrite Arm	Dendrite Arm	Dendrite Arm	Coating	Coating	Coating	Coating	Coating	
Company B Project 1 Project 2	3%													
	Coating	Coating	Coating	Coating	Coating	Particle Size	Particle Size	Particle Size	Particle Size	Particle Size	Particle Size	Particle Size	Particle Size	

The PAXCam Camera interface also received updates in the form of a cleaner, more organized, intuitive user-interface.

The foundation that made PAX-it such a successful tool still remains.

Electronic Filing System

- Simple organizational structure with easy navigation
- User may title and group folders
- Tools for security at different user levels
- Complete project management within project folders
- Image database searches are extensive and can go across all folders

◎·云·四·石·母·	N 2 9
a Select Fag Tile Cuttores	Mag.
All holisty	12 Fort Captor
ALL ROBINS	10/PeerCoupler
And and	10/free Cagle
and Nucleary	10:Per Crate
And And	10/Ren Cagle
Augusta .	10 Post Cuple
Robber	10/free Crupter
25 Maddate	10/Ren Cagler
Robber	10/Fort Cagle
	10 Peer Caster

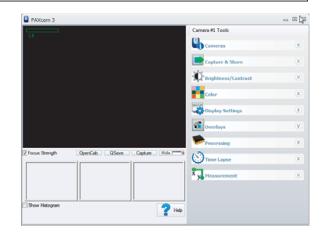
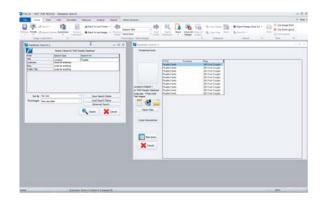




Image Database

- Search and retrieve data in Basic or Advanced modes
- Database fields that can be tailored to the user's needs
- Images, reports, video clips, data—any digital project file can be managed by PAX-it
- Search results show data table and image thumbnails
- ODBC-compliant database structure




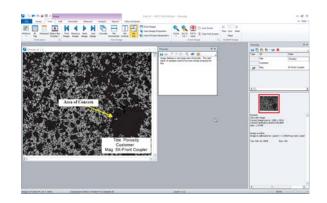


Image Processing

- Image stitching
- Image fusion—fuses multiple images of different focal planes into one
- Create video clips/AVI files
- Add editable annotations, voice narration, notes, and overlay templates


Measurement and Analysis

- New cost-effective modules available, streamlined to meet the customer's needs (reference 203-988-007)
- · From simple linear and area measurements to complex image analysis
- Metrology/Measurements on live preview of PAXCam (using optional Live Image Analysis Module)
- · Auto-thresholding of images based on density, color/grayscale, size, or shape
- Measurement on captured images

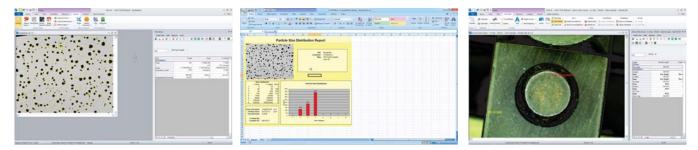

Networks and Communications

- Direct Email from within PAX-it
- Network licenses for all PAX-it functions and modules
- PAXnet for web-based sharing, downloading, and discussion
- Telemicroscopy modules for sharing remote views in real-time

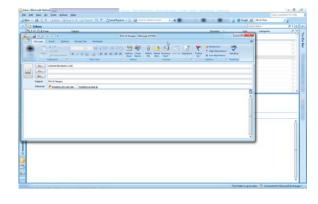

Specifications and part numbers may change. Consult LECO for latest information.

PAX-it™ and PAXcam™ are trademarks of Midwest Information Systems, Inc.

LECO Corporation 3000 Lakeview Avenue • St. Joseph, MI 49085 • Phone: 800-292-6141 • Fax: 269-982-8977 info@leco.com • www.leco.com • ISO-9001:2008 HQ-Q-994 • LECO is a registered trademark of LECO Corporation.



© 2015 LECO Corporation